



High bisulfite conversion efficiency using EpiMelt Bisulfite Modification Standard Kit on genomic DNA

Ea Marie Givskov Sørensen, Tina Ellegaard Kjeldsen, Lise Lotte Hansen

MethylDetect ApS, Aalborg, Denmark

Abstract

DNA methylation biomarkers are becoming increasingly relevant in assessing disease predisposition, diagnosis and detection of early-stage disease, personalized treatment, and post-treatment surveillance. In research, methylation analyses are often performed on limited or degraded biological material. Therefore, the sensitivity of the detection method is essential.

EpiMelt Methylation Detection Assays are based on the Methylation-Sensitive High-Resolution Melting (MS-HRM) technology, where the unique primer design and specific annealing temperature of the assay ensure robust and highly sensitive DNA methylation analysis.

In this application note, we show the performance of five EpiMelt Methylation Detection Assays and the EpiMelt Bisulfite Modification Standard Kit. We demonstrate this to be an efficient and sensitive method of bisulfite modification and methylation detection.

Introduction

DNA methylation is an epigenetic modification, which primarily consists of a methyl moiety attached to cytosines in CpG dinucleotides. In healthy cells, CpGs in gene-coding and non-gene-coding regions are generally methylated. At the same time, first exons, promoters, and enhancers containing high levels of CpG-content (CpG islands) typically are unmethylated^{1,2}. Hypomethylation is often a sign of active transcription, while hypermethylation is associated with gene transcription silencing³. However, the established DNA methylation patterns are disrupted in cancer, and CpG island methylation status at specific genes can have distinct consequences for patients at all stages of the disease.

One of the most common methods of investigating DNA methylation status is bisulfite conversion followed by a method of methylation detection. The bisulfite treatment converts unmethylated cytosines into uracils, while methylated cytosines are protected¹. Bisulfite conversion involves the use of harsh chemicals, which induces degradation of the DNA that then risks being lost during the subsequent DNA purification^{4,6}. However, the EpiMelt Bisulfite modification Standard Kit employs a bisulfite conversion step with high temperature and bisulfite concentration, and decreased processing time, a technique that efficiently and homogeneously converts DNA^{6,7}.

The EpiMelt Methylation Detection Assays are based on Methylation-Sensitive High-Resolution Melting (MS-HRM). MS-HRM is the methylation-independent amplification of the target region, followed by HRM⁸. During HRM, the fluorescence signal from a dsDNA intercalating dye is measured across an increasing temperature gradient. The melting profiles of the amplicons reflect whether the template DNA was methylated, as it will have a high GC content and therefore melts at a higher temperature, compared to amplicons from unmethylated template DNA⁹. The primers are designed to match the methylated template, to overcome PCR bias⁸. The EpiMelt Methylation Detection Assays contain a locus-specific primer mix and three ready-to-use controls (Methylation Positive Control, Assay Calibration Control, and Methylation Negative Control). The method is semiquantitative, and the methylation status of an unknown sample can be determined by comparison to the melting profiles of the three controls¹⁰.

In this application note, we demonstrate the robust and efficient bisulfite conversion and clean-up of formalin-fixed paraffin-embedded (FFPE) non-small-cell lung cancer samples, and the sensitive methylation detection using the EpiMelt Bisulfite modification Standard Kit in combination with the EpiMelt methylation detection assays targeting *BRCA1*, *BRCA2*, *HORMAD1*, *MGMT1*, and *MLH1*.

Materials

Reagents and consumables

Genomic DNA (gDNA) DNA was extracted from 5 × 10 μm from non-small-cell lung cancer (NSCLC) formalin-fixed paraffin-embedded (FFPE) sections using the High Pure FFPE DNA Isolation Kit (Roche, Germany) according to the manufacturer's protocol and stored in TE buffer. Samples were divided into low, medium, and high-quality DNA based on an agarose gel electrophoresis.

Materials

Qubit™ 1 × dsDNA HS Assay Kit (Thermo Fisher Scientific, cat. #Q33230).
 EpiMelt Bisulfite Modification Standard Kit (MethylDetect ApS, cat. # EPI-BIS-200).
 EpiMelt *BRCA1* test assay (MethylDetect ApS).
 EpiMelt *BRCA2* test assay (MethylDetect ApS).
 EpiMelt *HORMAD1* test assay (MethylDetect ApS).
 EpiMelt *MGMT* test assay (MethylDetect ApS).
 EpiMelt *MLH1* test assay (MethylDetect ApS).
 EpiMelt Real-Time PCR Master Mix (MethylDetect ApS, cat. # EPI-qPCR-200).
 LightCycler® 480 Multiwell plate 96, white (Roche, cat. # 04729692001).

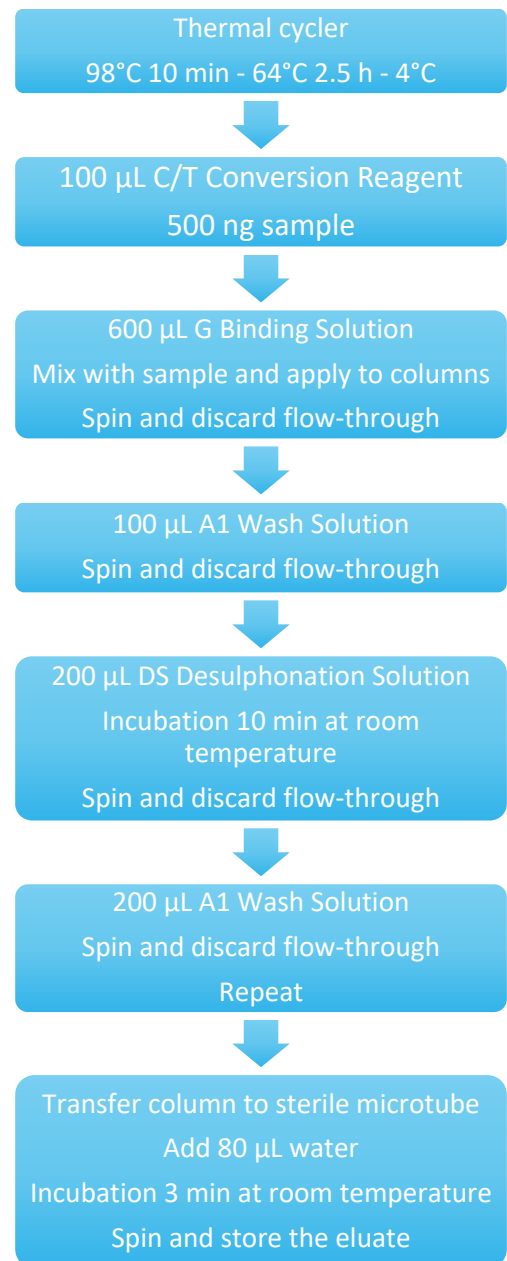


Figure 1: Bisulfite conversion and purification workflow.

Instruments

Qubit™ 3 Fluorometer (Thermo Fisher Scientific, cat. # Q33216).

LightCycler® 480 System (Roche, cat. # 05015278001).

Methods

Bisulfite conversion and purification

Prior to bisulfite conversion and DNA purification, the gDNA concentration was measured using the Qubit™ 1 × dsDNA HS Assay on the Qubit™ 3 Fluorometer.

Water and D Dilution Buffer were added to the C/T Conversion Reagent from the EpiMelt Bisulfite Modification Standard Kit to solubilize the solid crystalline. For bisulfite conversion, 500 ng gDNA from each of the three samples was diluted with water to a volume of 50 µL and used as input. Bisulfite treatment and DNA purification were performed using the EpiMelt Bisulfite Modification Standard Kit according to the workflow depicted in Figure 1.

Amplification and high-resolution melting

Five EpiMelt Methylation Detection Assays were used to analyze the gDNA samples, targeting *BRCA1*, *BRCA2*, *HORMAD1*, *MGMT*, and *MLH1*. Each EpiMelt Methylation Detection Assay contains a locus-specific primer mix and three ready-to-use controls. The EpiMelt Controls or the converted and purified gDNA samples were combined with 100 µL EpiMelt Real-Time PCR Master Mix, 3.0 µL water, and 1.0 µL EpiMelt Primer Mix according to Table 1.

Reagent	Volume
EpiMelt Real-Time PCR Master Mix 2 ×	10.0 µL
Nuclease-free water	3.0 µL
EpiMelt Primer Mix (10 mM)	1.0 µL
EpiMelt Control (5ng/mL) or converted, purified gDNA sample	6.0 µL
Total	20.0 µL

Table 1: PCR mix.

The PCR amplification and subsequent high-resolution melting were performed on the LightCycler® 480 Instrument II in LightCycler® 96-well plates using the setting Detection format » SYBR Green I/HRM dye».

PCR program	Cycles	Temperature (°C)	Acquisition mode	Hold (sec)	Ramp rate (°C/sec)	Acquisitions (per °C)
Pre-incubation	1	95	None	600	04.04	-
Amplification	50	95	None	15	04.04	-
		<i>BRCA1</i> : 56 <i>BRCA2</i> : 59 <i>HORMAD1</i> : 50 <i>MGMT</i> : 59 <i>MLH1</i> : 60	None	10	02.02	-
		72	Single	15	04.04	-
High-resolution melting	-	95	None	15	04.04	-
		60	None	60	02.02	-
		95	Continuous	-	00.01	50

Table 2: PCR and HRM program. The annealing temperature is specific to the assay and may vary between platforms.

The LightCycler® 480 Instrument II program for PCR amplification and the high-resolution melting are shown in Table 2. The listed annealing temperature is dependent on the assay. An annealing temperature of 56°C was used for EpiMelt *BRCA1*, 59°C for EpiMelt *BRCA2*, 50°C for EpiMelt *HORMAD1*, 59°C for EpiMelt *MGMT*, and 60°C for EpiMelt *MLH1*. No template controls were included for each assay, and all reactions were analyzed in triplicates. Data were analyzed using the LightCycler® 480 Software.

Results

The use of the EpiMelt Methylation Detection Assays and the EpiMelt Bisulfite Modification Standard Kit resulted in efficient bisulfite modification of all nine samples, followed by sensitive methylation detection. This is illustrated by the amplification curves (Figure 2A - E) and the relative signal difference (-d/dT) plots (Figure 2F - J) for each assay, and the Ct values for each sample and the EpiMelt Controls (Table 3).

The sample amplification curves were steep and reproducible for all EpiMelt Methylation Detection Assay replicates (Figure 2A-I). The high amplification signals and low Ct values (Table 3) suggest excellent template recovery and preservation of DNA integrity. In the relative difference plots, the peaks are uniform, and all three controls are clearly separable (Figure 2F-J). The melting profiles of the samples suggest complete bisulfite conversion, as they display uniform melting profiles that are easily comparable to the controls.

Figure 2.

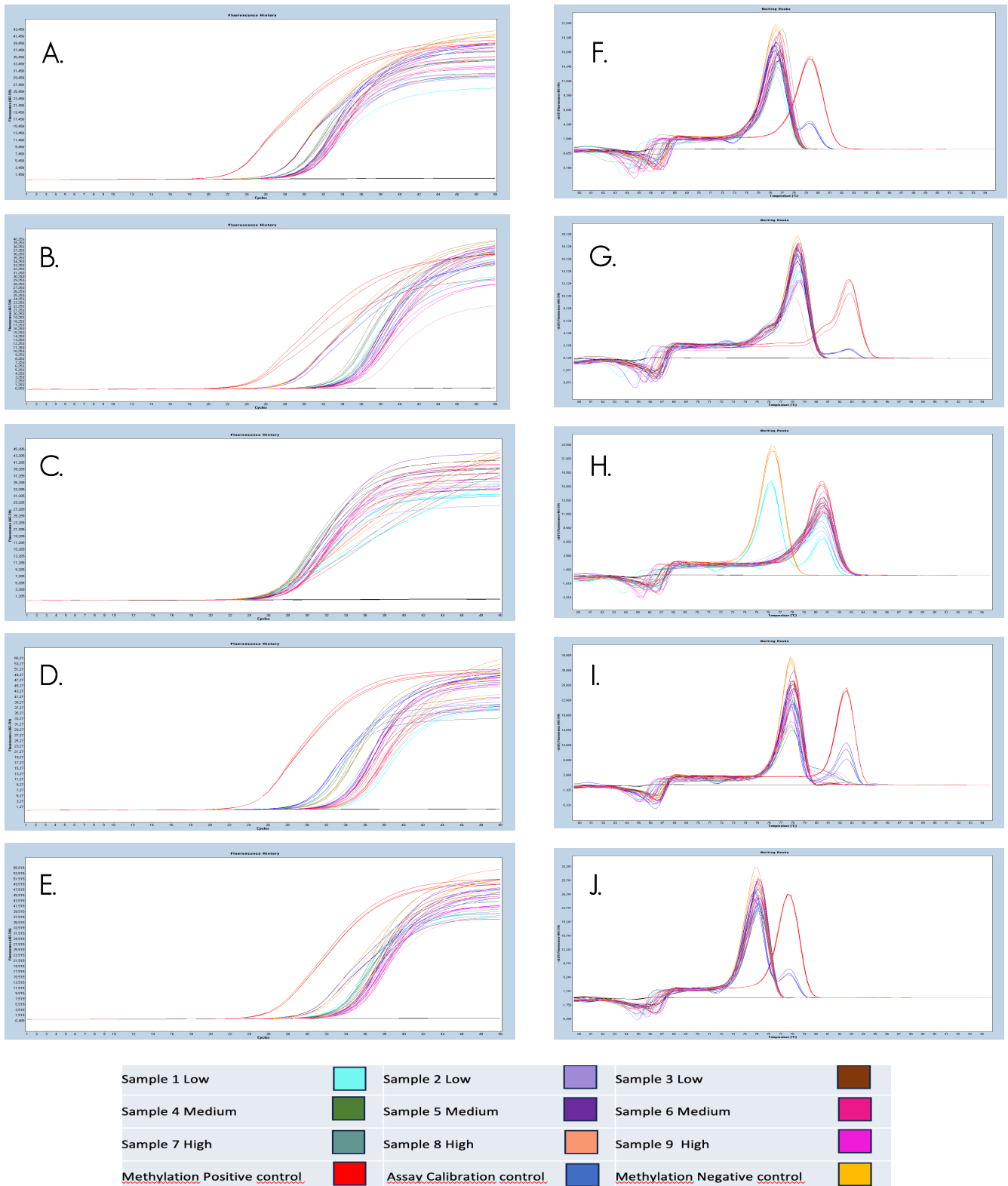


Figure 2: Amplification of nine FFPE DNA NSCLC samples and EpiMelt Controls using the EpiMelt *BRCA1* Assay (A), EpiMelt *BRCA2* Assay (B), EpiMelt *HORMAD1* Assay (C), EpiMelt *MGM1* Assay (D), and EpiMelt *MLH1* Assay (E). Melting profiles of nine FFPE DNA

NSCLC samples and EpiMelt Control PCR products from EpiMelt *BRCA1* Assay (F), EpiMelt *BRCA2* Assay (G), EpiMelt *HORMAD1* Assay (H), EpiMelt *MGMT* Assay (I), and EpiMelt *MLH1* Assay (J).

	<i>BRCA1</i>	<i>BRCA2</i>	<i>HORMAD1</i>	<i>MLH1</i>	<i>MGMT</i>
Methylation Positive Control	22	24	27	25	24
Assay Calibration Control	26	27	26	31	29
Methylation Negative Control	26	27	26	31	34
Sample 1 Low Quality	30	34	28	33	35
Sample 2 Low Quality	30	34	28	34	33
Sample 3 Low Quality	30	34	28	34	34
Sample 4 Medium Quality	29	34	27	33	32
Sample 5 Medium Quality	30	35	28	34	34
Sample 6 Medium Quality	31	35	28	34	34
Sample 7 High Quality	29	34	27	32	30
Sample 8 High Quality	30	36	28	34	34
Sample 9 High Quality	30	35	28	34	34

Table 3: C_t values for the EpiMelt Controls and each gDNA sample for the five EpiMelt Methylation Detection Assays targeting *BRCA1*, *BRCA2*, *HORMAD1*, *MLH1*, and *MGMT*.

Conclusion

In this application note, we demonstrate that the use of the EpiMelt Bisulfite Modification Standard Kit leads to efficient and complete bisulfite conversion of genomic DNA extracted from blood samples, with excellent template recovery. This is obtained with support from the EpiMelt Methylation Detection Assays, showing state-of-the-art methylation sensitivity.

References

- 1 Clark, S. J., Statham, A., Stirzaker, C., Molloy, P. L. & Frommer, M. DNA methylation: Bisulphite modification and analysis. *Nature protocols* **1**, 2353-2364 (2006). <https://doi.org/10.1038/nprot.2006.324>
- 2 Nishiyama, A. & Nakanishi, M. Navigating the DNA methylation landscape of cancer. *Trends Genet* **37**, 1012-1027 (2021). <https://doi.org/10.1016/j.tig.2021.05.002>
- 3 Unnikrishnan, A. *et al.* The role of DNA methylation in epigenetics of aging. *Pharmacol Ther* **195**, 172-185 (2019). <https://doi.org/10.1016/j.pharmthera.2018.11.001>
- 4 Miura, F., Enomoto, Y., Dairiki, R. & Ito, T. Amplification-free whole-genome bisulfite sequencing by post-bisulfite adaptor tagging. *Nucleic Acids Research* **40**, e136-e136 (2012). <https://doi.org/10.1093/nar/gks454>
- 5 Tanaka, K. & Okamoto, A. Degradation of DNA by bisulfite treatment. *Bioorganic & Medicinal Chemistry Letters* **17**, 1912-1915 (2007). <https://doi.org/10.1016/j.bmcl.2007.01.040>
- 6 Yi, S., Long, F., Cheng, J. & Huang, D. An optimized rapid bisulfite conversion method with high recovery of cell-free DNA. *BMC Molecular Biology* **18**, 24 (2017). <https://doi.org/10.1186/s12867-017-0101-4>
- 7 Genereux, D. P., Johnson, W. C., Burden, A. F., Stöger, R. & Laird, C. D. Errors in the bisulfite conversion of DNA: modulating inappropriate- and failed-conversion frequencies. *Nucleic Acids Res* **36**, e150 (2008). <https://doi.org/10.1093/nar/gkn691>
- 8 Wojdacz, T. K., Hansen, L. L. & Dobrovic, A. A new approach to primer design for the control of PCR bias in methylation studies. *BMC Res Notes* **1**, 54 (2008). <https://doi.org/10.1186/1756-0500-1-54>
- 9 Hussmann, D. & Hansen, L. L. Methylation-Sensitive High Resolution Melting (MS-HRM). *Methods in molecular biology (Clifton, NJ)* **1708**, 551-571 (2018). https://doi.org/10.1007/978-1-4939-7481-8_28
- 10 Wojdacz, T. K. & Dobrovic, A. Melting curve assays for DNA methylation analysis. *Methods in molecular biology (Clifton, NJ)* **507**, 229-240 (2009). https://doi.org/10.1007/978-1-59745-522-0_17